

# Turn in-house FR-X OFF

---

## FR-X

Decrease gradually the voltage and intensity from 45 kV and 66 mA to 20 kV and 10 mA. First step 66→60 mA, then 45→40 kV. Wait 10 s. Then decrease by steps of 10. (Lower the intensity first)

Turn X-RAY OFF on the XG Control panel. Wait 1h.

Stop the Vacuum on the XG Control panel. Wait 1h.

POWER OFF on the FR-X main panel.

Switch OFF at the back of the FR-X.

Switch the main power source OFF.

---

## Cryostream

Program / Ramp (change with dial) / Enter

Choose the speed (up to 360K/h) / Enter Target / Ambient / Enter Enter / Enter Start

When ambient temperature is reached, stop Turn OFF at the back

---

## Engine Room

Switch OFF the nitrogen compressor in the engine room using the touch panel.

Switch OFF the main switch of the other big device in the engine room.

From:

<https://bsi.inscog.eu/> - **BSI wiki**

Permanent link:

<https://bsi.inscog.eu/doku.php?id=crystallography:collection:frx:off&rev=1469103994>

Last update: **2023/11/01 20:17**

