

# Setup an automated data collection at Glacios

## Beam microprobe and nanoprobe calibrations

- Introduce the Cross Grating Grid in the column.
- Go to x34 Imaging status.
- Select 2 positions: a carbon square and an empty square.
- Introduce the Condenser 2 of 30 um
- Go to x280 Imaging status.
- In DigitalMicrograph click View and place the mouse arrow on a feature.
- Use Alpha Toggle 20° in Stage<sup>2</sup> window (TEM) to rock the grid.
- Correct with **Z axis knobs** on the console.
- Repeat a second time.
- Go to x45k Imaging status
- In Low Dose Control window (SEM) click on Go to: Record.
- Push the **Eucentric Focus** knob on the console.
- Inset the screen (FC).
- Push the **Wobbler** knob on the console.
- Adjust the Z height on the screen with the **Z axis knobs** on the console.

## Microprobe

Focus the beam and center it (with the trackball).

- Go to microprobe (TEM)
- Push the **Eucentric Focus** knob on the console.
- In the **Direct Alignments** window:

1. Adjust the [Beam tilt pp X](#)
  - Click on **Beam tilt pp X**
  - Adjust with **Multifunction X/Y** knobs
  - Click **Done** at the bottom of the window.
2. Adjust the [Beam tilt pp Y](#)
  - Click on **Beam tilt pp Y**.
  - Adjust with **Multifunction X/Y** knobs.
  - Click **Done** at the bottom of the window.
3. Adjust the [Beam shift](#)
  - Click on **Beam shift**.
  - Adjust with **Multifunction X/Y** knobs.
  - Click **Done** at the bottom of the window.
4. Adjust the [Rotation center](#)

- Click on **Rotation center**.
- Adjust with **Multifunction X/Y** knobs.
- Click **Done** at the bottom of the window.

## Nanoprobe

Focus the beam and center it (with the trackball).

- Go to nanoprobe (TEM)
- Push the **Eucentric Focus** knob on the console.
- Check the Condenser 2 centering and adjust if necessary.
- In the **Direct Alignments** window:
  1. Adjust the [Beam tilt pp X](#)
    - Click on **Beam tilt pp X**
    - Adjust with **Multifunction X/Y** knobs
    - Click **Done** at the bottom of the window
  2. Adjust the [Beam tilt pp Y](#)
    - Click on **Beam tilt pp Y**
    - Adjust with **Multifunction X/Y** knobs.
    - Click **Done** at the bottom of the window
  3. Adjust the [Beam shift](#)
    - Click on **Beam shift**
    - Adjust with **Multifunction X/Y** knobs
    - Click **Done** at the bottom of the window
  4. Adjust the [Rotation center](#)
    - Click on **Rotation center**
    - Adjust with **Multifunction X/Y** knobs
    - Click **Done** at the bottom of the window

[After all calibrations save the settings](#)

- Menu: *Settings* > *Save*
- In **Low Dose Control** window check **Continuous update**
- Click on Update it
- Uncheck **Continuous update**
- Menu: *Settings* > *Save*

## Correct astigmatism

- Verify that the **Turbo Auto Off** Vacuum option is checked (TEM)

- Go to x45k Imaging status
- In the Camera and script window (SEM) open the **Setup**
- In Record Tab, set 2 sec exposure, processing unnormalized, align frames, save frames and the objective diaphragm out
- If the signal is too weak, increase the time, concentrate the beam or change the carbon area
- **Correct astigmatism by CTF**
  - Menu: Calibration > Focus & Tuning > Correct astigmatism by CTF
  - Open FFT window with Shift+F
  - Repeat several times until the corrected parameters have converged
- **Coma free align by CTF**
  - Menu: Calibration > Focus & Tuning > Coma free align by CTF
  - Open FFT window with Shift+F
- **Coma vs image shift**
  - Menu: Calibration > Focus & Tuning > Coma vs image shift
  - Open FFT window with Shift+F
  - If the signal is too weak, increase the time, set binning, concentrate the beam or change the carbon area
  - parameters for R 1.2/1.3 grids: 5  $\mu\text{m}$
  - parameters for R 2/2 grids: 6  $\mu\text{m}$

From:

<https://bsi.inscog.eu/> - **BSI wiki**

Permanent link:

[https://bsi.inscog.eu/doku.php?id=glacios\\_collect&rev=1644592817](https://bsi.inscog.eu/doku.php?id=glacios_collect&rev=1644592817)

Last update: **2023/11/01 20:14**

