


Make sure Microscope is in good shape (ppt associated)

- a. Make sure HT and FEG are on
 - On TEM User Interface (TUI)
 - setup tab > High Tension panel High tension button is highlighted yellow
 - setup tab > FEG Control panel Operate button is highlighted yellow

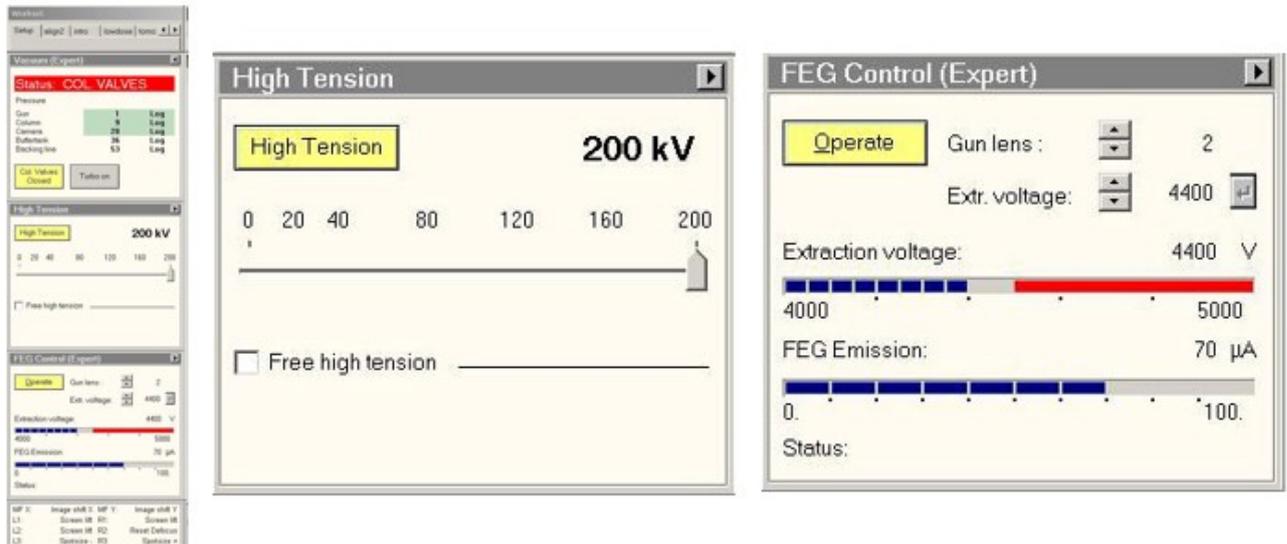

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
- b. Make sure column valves are closed and turbo pump is on
 - On TEM User Interface (TUI)
 - setup tab > Vacuum panel Col. Valves closed button is highlighted yellow
 - setup tab > Vacuum panel Turbo Pump button is highlighted yellow

IMAGE

- c. Make sure that EM column vacuum is stable (IGP1 = 6)
- setup tab > Vacuum panel Column value must be 6 IMAGE

- Setup the microscope for recording the Atlases

- On TEM user interface (**TUI**).



From:

<https://bsi.inscog.eu/> - **BSI wiki**

Permanent link:

https://bsi.inscog.eu/doku.php?id=tf20_screening&rev=1674236757

Last update: **2023/11/01 20:16**

